Aseptic Sampling Valves

Sanitary
flow
equipment

Tipo **VPAK**

Pres campion asettica

CARATTERISTICHE

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
STERILE SAMPLING

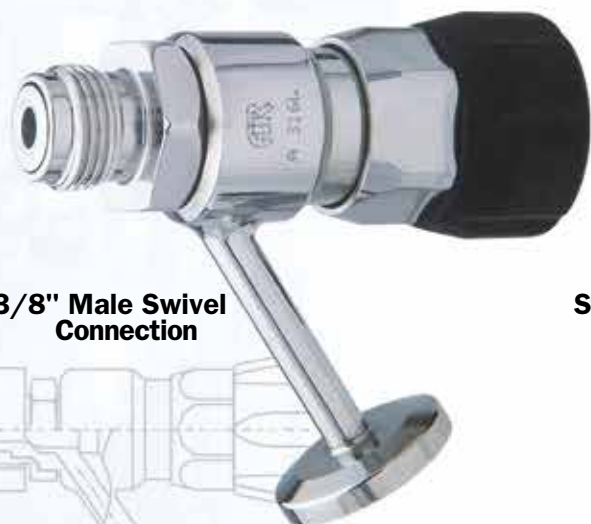Developed for taking samples in a sterile way, the special configuration assures effective cleaning (CIP) and sterilization (with saturated steam) of product contact surfaces of valve and seating plug with the valve in closed position. Once installed, the sealing plug shuts off flush with the inlet port connection with a zero dead leg.

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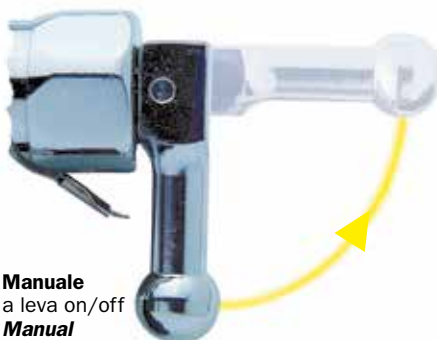
Manuale
a leva on/off
Manual
stainless steel on-off lever

Pneumatico inox on/off
Air operated stainless steel air-spring actuators

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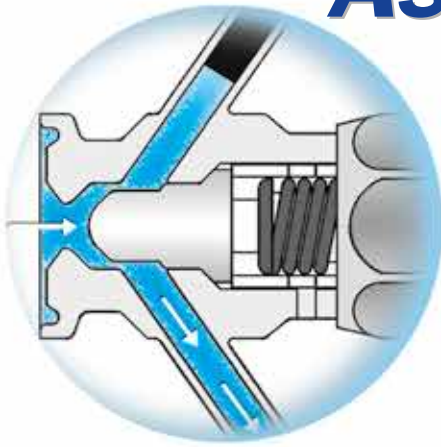
DOCUMENTATION

All valves are fully traceable for validation processes. Each valve is accompanied by material test certificates EN 10 204 3.1B 3A standards compliance certificates, FDA regulations compliance certificates, and tapes detailing the valve interior surfaces Ra.

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
IMPORTANT!

For reliable performance there must be a safety margin between the spring load of the seating plug and the maximum pressure in the product line. Higher pressures will force open the valve. Please indicate the maximum pressure of your applications.

Sampling phase

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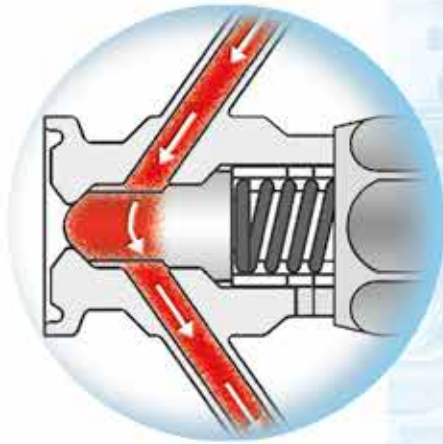
fig. 1
Valvola aperta (fase di prelievo)
Valve open (sampling phase)

SIP phase

■ Vapore
Steam

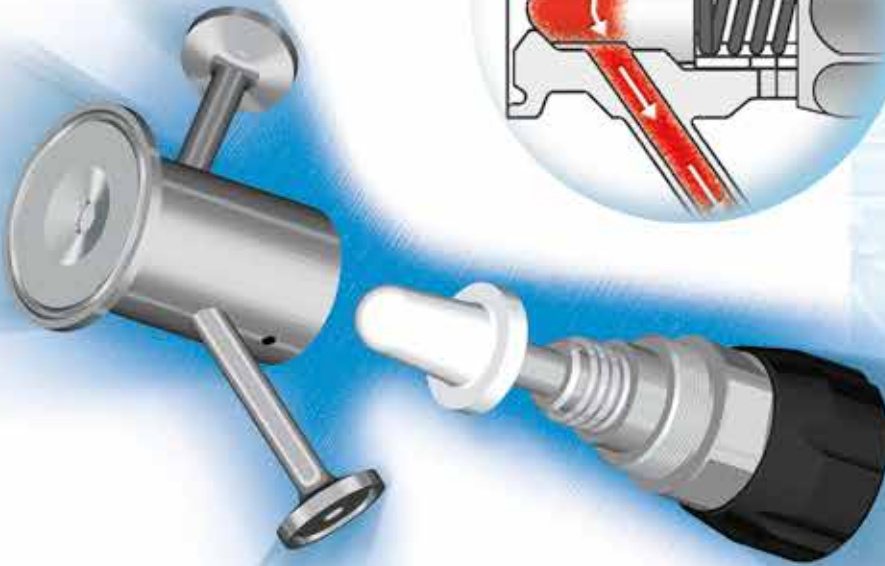fig. 2
Valvola chiusa (fase di CIP/steriliz.)
Valve closed (CIP/steril. phase)

VPA-VPAK Ordering information

To specify the part completely, start with the product description and select the additional options as shown below :

	<i>Actuator type</i>		<i>Sample port</i>		<i>diaphragm</i>	<i>Surface finish</i>				
	manual knob M		clamp UTC		silicone X	Ra 0,5 micron as standard				
	manual lever ML		butt weld USD		TFM 1600 PTFE G	EP electro-polished on request				
	pneumatically actuated P		hose barb UPG							
	pneumatic & switch PS		Staubli conn. UST							
<i>Valve type</i>	VPAK	M	ETC	UTC	VTC	G	1.4404	Ra		Atex Ex II 2 GD T4
	VPA									
			<i>Process connection</i>		<i>CIP/SIP port</i>	<i>Material</i>				
			clamp ETC		VTC clamp VTC	1.4404 as standard				
			butt weld ESD		VSD butt weld VSD	1.4435 on request				
			Ingold connection EIN		VPG hose barb VPG					
			BSP thread EGM							
			flanged EFL							

Aseptic sampling valve

Type **VPA**

CHARACTERISTICS

The **AR** sampling valve range has been designed for sterile sampling applications in the Chemical and Pharmaceutical industries. Its compact design allows easy installation on small vessels or process lines.

FULLY AUTOCLAVABLE,

Valve body and head are made out of certified 316L bar stock.

Damage to the seating plug and unexpeted material in the outlet are indications of overpressure. A leak detector hole is placed immediately behind the outlet.

On request the valve can be equipped with a CIP system.

SURFACE FINISH

External mechanical "mirror" polish
Internal mechanical polish Ra < 0.4 µm

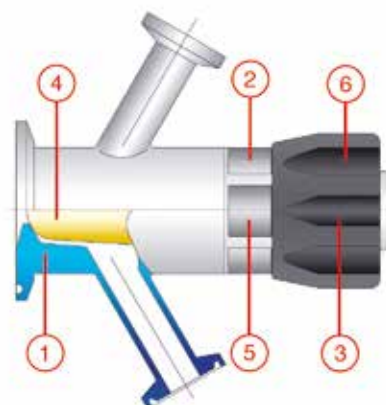
VALVE BODY connections available in:

- weld ends
- clamp ends

Other types available upon request.

MATERIALS

- 1 AISI 316L valve body
- 2 AISI 316L valve head
- 3 AISI 316L stem
- 4 Silicone or PTFE membrane
- 5 AISI 304 spring
- 6 adjustable autoclavable PTFE knob

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Pressione: 8 bar
Temperatura: -30°C a +150°C
Portata: 1.200 lt/h a 2 bar
Membrana: silicone - PTFE

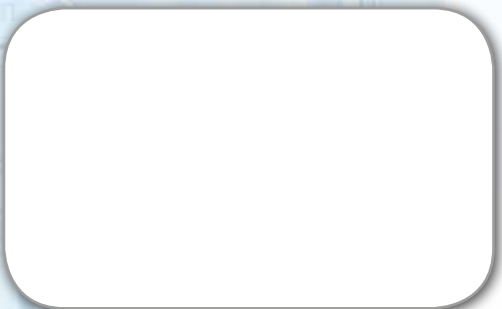
MAX WORKING CONDITIONS

Pressure: 8 bar (116 Psi)
Temperature: -30°C up to +150°C
(-22°F up to +302°F)
Flow rate: 1.200 lt/h at 2 bar
(317 US gal/hr at 29 Psi)
Diaphragm: silicone - PTFE

Sanitary flow equipment

AERRE INOX s.r.l.
Via Gerola, 4
I-26010 Fiesco (Cremona)
Tel. +39 0374 370 828 Fax +39 0374 370 833
www.aerreinox.it

Get the information you need and more at info@erreinox.it