# Sanitary Sampling Valves

Sanitary  
flow  
equipment

# Rps Available Body

**Standard  
CLAMP 1/2" BS**

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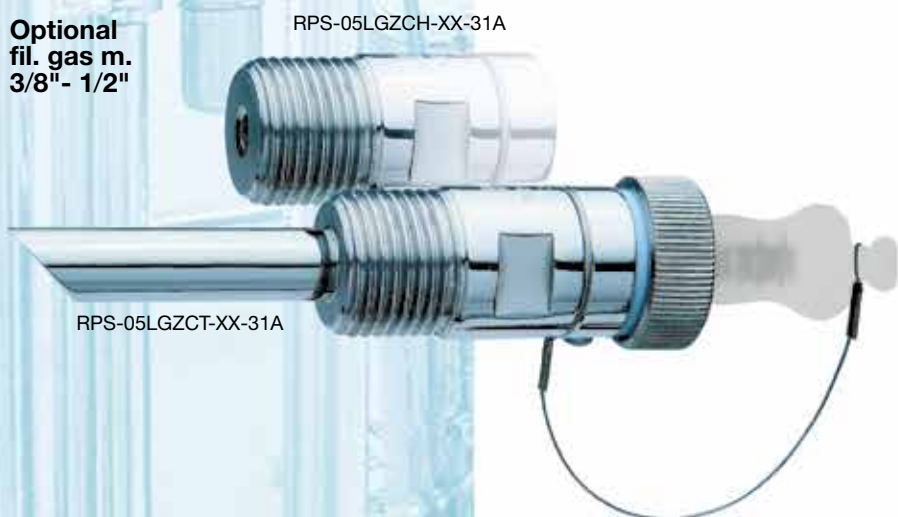
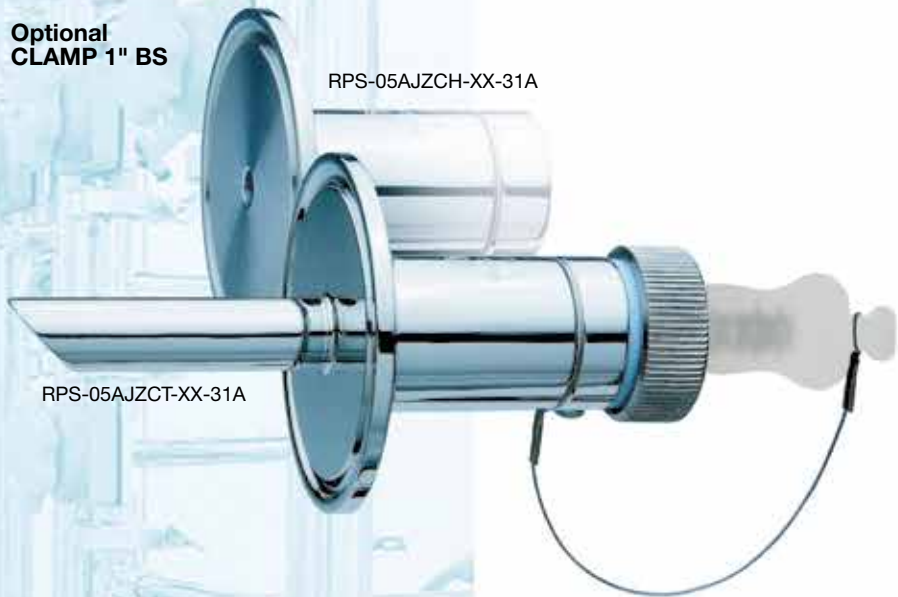
RPV-05AGZCH-XX-31A


RPV-05AGZGH-XX-31A

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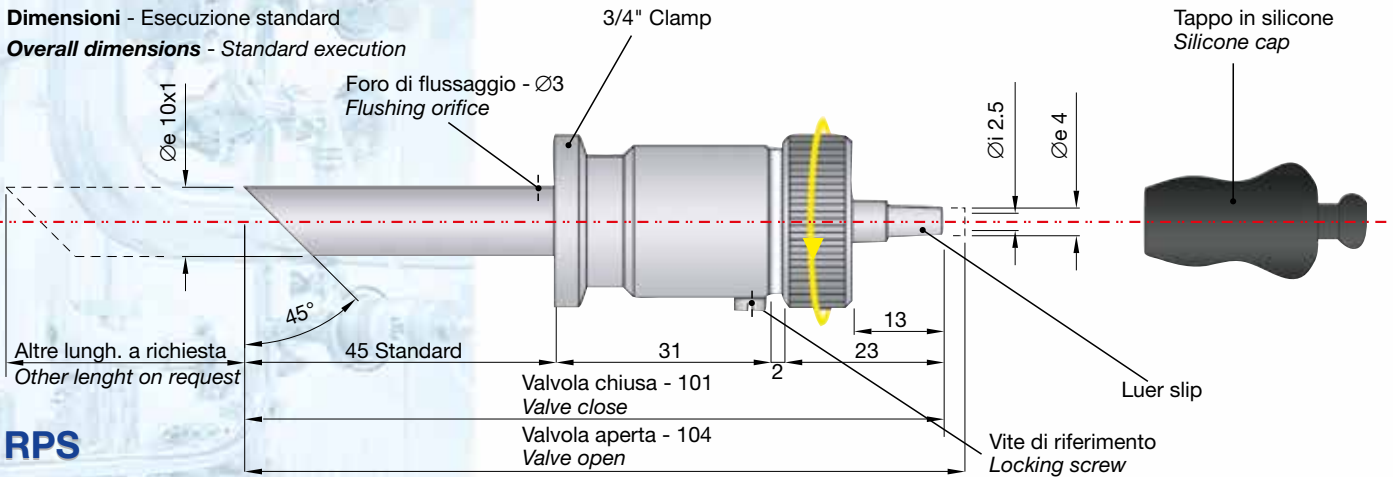
Overall dimensions - Standard execution

RPS

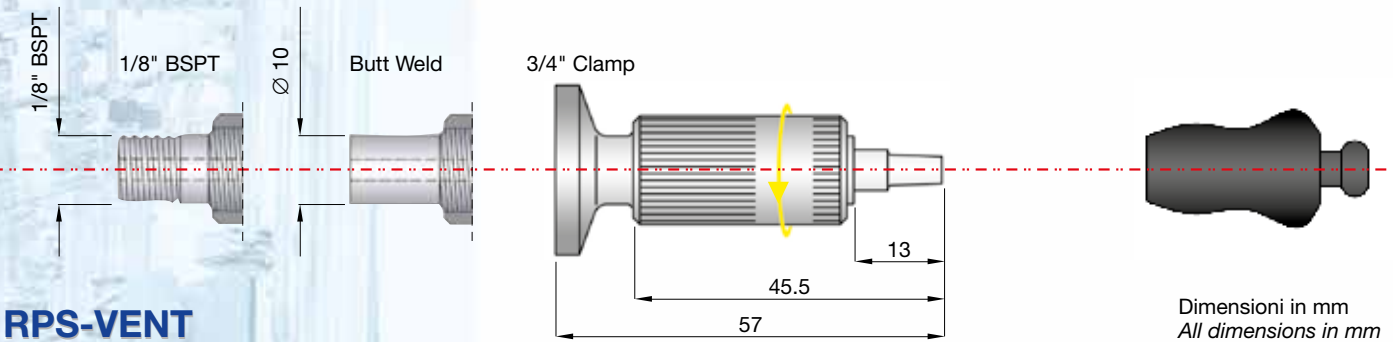RPS-VENT

Dimensioni in mm  
All dimensions in mm

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## INSTALLATION

Depending on the type of connection, the valve may be installed by welding, screwing or clamping the body.  
As the direction of flow is one-way, check the direction of the reference screws to ensure the body valve flow is correct.  
For all types of connection, ensure the luer outlet is horizontal to the valve and parallel with the floor.  
If flow is vertically upwards, the reference screws must be placed at the bottom, while if the flow is vertically downwards, the reference screws must be placed at the top.

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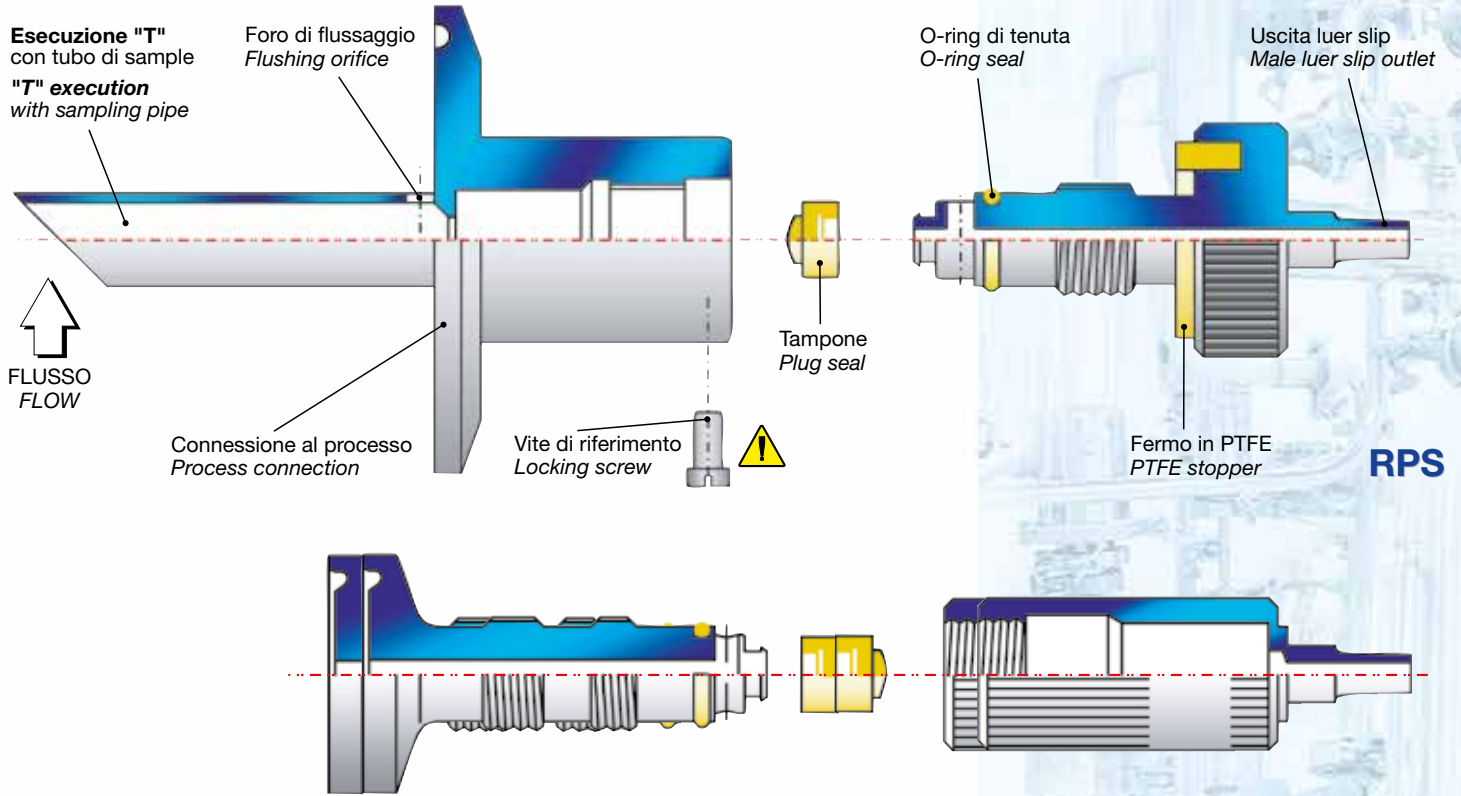
## DOCUMENTATION

All valves are fully traceable for validation processes.  
Each valves is accompanied by material test certificates EN 10 204 3.1 cGMP standards compliance certificates, FDA regulations compliance certificates, and tapes detailing the valve interior surfaces Ra.

# Sanitary sampling valve

## Ricambi Spare parts

### cod. KPS-05-VV

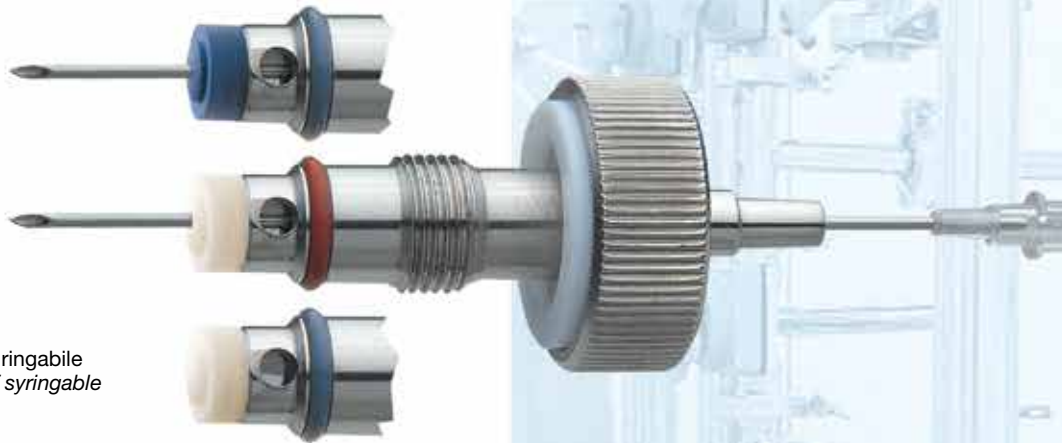
Tamponi e o-ring in FKM - siringabile  
Plug seal and o-ring in FKM - syringable

### cod. KPS-05-XX

Tamponi e o-ring in SILICONE - siringabile  
Plug seal and o-ring in SILICONE - syringable

### cod. KPS-05-TV

Tamponi in TFM PTFE e o-ring in FKM - NON siringabile  
Plug seal in TFM PTFE and o-ring in FKM - NOT syringable

## RPS Ordering information

To specify the part completely, start with the product description and select the additional options as shown below:

	<b>AG</b>	1/2" clamp						
	<b>AJ</b>	1" clamp						
	<b>QC</b>	butt weld						
<b>RPS</b>	<b>LC</b>	1/4" BSP						
<b>RPQ</b>	<b>LE</b>	3/8" BSP	<b>H</b>	flat face				
<b>RPV</b>	<b>LG</b>	1/2" BSP	<b>T</b>	deep tube				
							surface standard finish	
								material AISI 316L 1.4404

<b>RPS 05 AG ZC H XX 31 A</b>	
HOLE	<b>ZC</b> luer slip
	<b>PA</b> rubber hose
	<b>ZG</b> quick connection
	<b>XX</b> silicone seals
	<b>VV</b> Fkm seals
	<b>TV</b> Tfm/Fkm seals

# Sanitary sampling valves

## CHARACTERISTICS

The **RPS** fully autoclavable sampling valves are designed for aseptic sampling of liquids and pharmaceutical products.

Made entirely in **AISI 316L**, the standard version is in the form of a rod with a 1/2" clamp and no welding.

On request, the connection to the process may be modified as indicated.

Easily installed onto distribution lines or onto tanks, this valve can be kept in aseptic conditions between one sample and another by using a silicon cap which closes the "luer slip" outlet. It's special design prevents deposits of bacteria and particles, possibly distorting the results of bacteriological tests carried out while allowing samples to be taken from the middle of the line enabling a representative sample to be taken. Extremely economical and very easy to use, the valve opens by turning the head knurled ratchet in an anticlockwise direction and then clockwise to return it to the closed position.

## TECHNICAL DATA

### MATERIALS

Valve body: **AISI 316L**

Shaft: **AISI 316L**

O-ring: Standard silicone  
Optional FKM, EPDM

Plug: Standard silicone  
Optional FKM or TFMPTE

### MAX WORKING CONDITIONS

Max Pressure: to 10 bar and 150°C

### PROCESS CONNECTIONS

Standard: clamp 1/2" BS

Optional: clamp 1" BS

3/8"-1/2" gas male thread  
butt-weld

### OUTLET CONNECTIONS

Standard: LUER SLIP joint

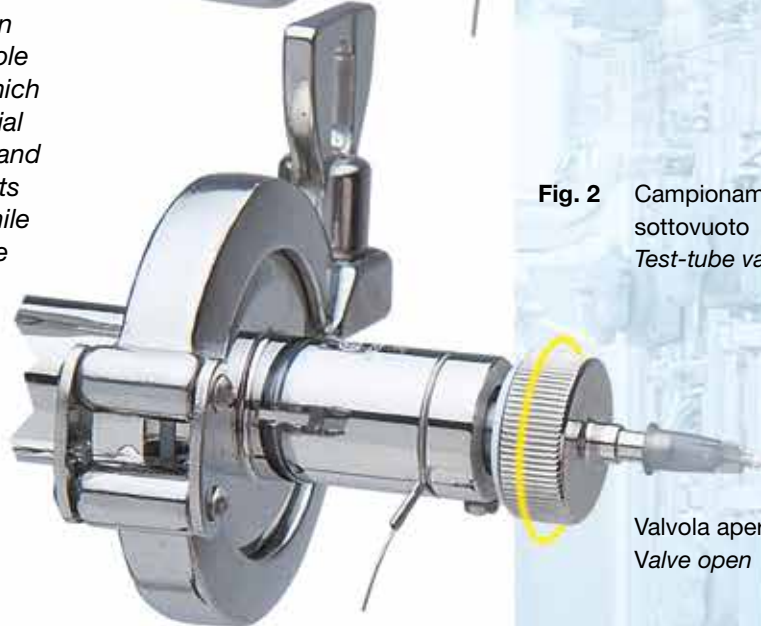
Optional: Rubber hose Ø 6mm  
Rapid connection

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